

Job Safety Analysis

Safety Information for The University of North Carolina at Chapel Hill

JOB SAFETY ANALYSIS TOPIC:

Sanitizing IVC Full Racks with the Rack Washer



	Title	Work Task	Hazards	Controls				
	Preplan the work	Gather proper tools and equipment	Improper tools					
		Gather PPE: Personal Protective Equipment (PPE) consisting of yellow	Improper/No Equipment Inspection					
		isolation gown, gloves, steel toe shoes	Improper PPE inspection/selection					
	Before the job	Don necessary PPE Inspect Rack Washer Door, Washer area and the IVC Full rack. Make sure Rack Washer is powered on and ready for service.	Improper Donning	Improperly donning or not donning PPE can lead to an unnecessary exposure. Familiarize yourself with the location and proper use of all controls and safety devices. Make sure the rack washer is free of obstructions such as: racks, equipment, cages etc.				

	 Open the rack washer door. Push the IVC Full Rack inside the Rack Washer. Close the washer door. Start Cycle. 		Ergonomic Fall Hazards Foot Hazard Hand injury	Check throughout the cycle to make sure the rack has not moved or adjusted during the cleaning cycle.		
Completing the sanitizing operation	Open the washer door. Remove the full rack. Doff PPE and dispose of it in proper locations		Ergonomic Fall Hazards Foot Hazard Hand injury	Clean up any excess water or spills immediately to prevent falling. Check the IVC rack for any damage after removing from the washer. (Report any damages to the supervisor and remove the rack from service use)		
Training	Employees must have hands on training in the proper use of this piece of equipment.		_			
Created By	Neah Tucker, Facilities Safety Officer Derick Pacio, Operations Manager					

Contact Info

For more information about this JSA and other JSAs, contact:

UNC-CH Department of Environment, Health and Safety

1120 Estes Drive Extension, Chapel Hill NC 27599 CB# 1650

(919) 962-5507 **#** https://ehs.unc.edu/



